750 x 1500mm, Low Profile Lab Table

#33-782

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General

Type:	Metric
Vibration Rating:	5 Hz: Vertical = 70 - 85%, Horizontal = 75 - 90% 10 Hz: Vertical = 90 - 97%, Horizontal = 90 - 97%


Physical & Mechanical Properties

Thickness (mm):	50.40
Length (mm):	1,500.00
Width (mm):	750.00
Construction:	Top: Stainless Steel Core: All Steel, High Density, Cell-Sized Honeycomb Core Frame: Medium Texture Black Powder Coat
Feed Pressure (psi):	80 (nitrogen or air)
Flatness (inches):	±0.005
Height (inches):	31.00
Load Capacity (lbs):	Gross: 1400 Net: 350
Weight (lbs):	600.00


Threading & Mounting

Mounting Threads:	M6 on 25 Centerline
-------------------	---------------------

Damping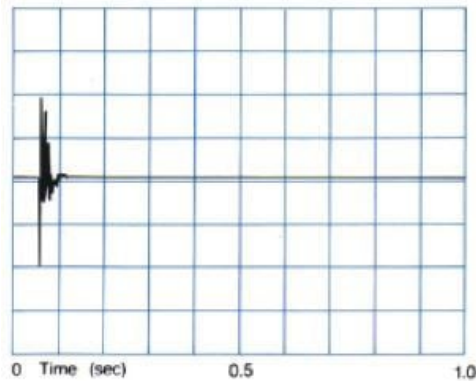

Compliance

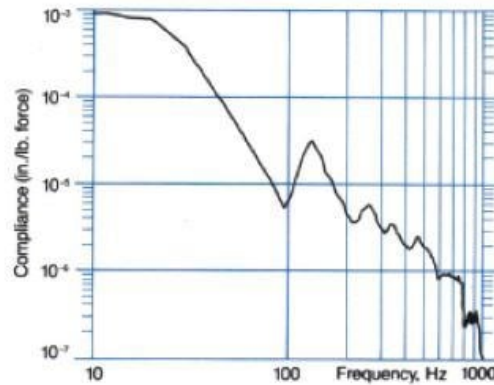
CleanBench™ Steel Top — 30 x 48 x 2 in. (750 x 1200 x 50 mm)

Structural Damping determines how quickly an excited resonance in a table top decays. The simplest way to measure damping is to hit the table top with a hammer and measure the decay with an accelerometer and oscilloscope or spectrum analyzer. The height of a resonance peak in the "compliance curve" also measures damping.

Compliance is a reciprocal measurement of the dynamic stiffness of a table top. The data are obtained by inputting a measured force to the table top with a calibrated hammer and measuring the resultant acceleration (or displacement) with an accelerometer. Compliance is the ratio of displacement to force expressed as a function of frequency.

Isolator Performance

All CleanBench tables incorporate TMC's Gimbal Piston Air Isolators as a standard feature. The Gimbal Piston has consistently shown to out-perform other air isolators in side-by-side testing. It offers outstanding low frequency vibration isolation in all axes and maintains its performance specifications even when subjected to extremely low input levels of excitation. Proprietary damping techniques allow the Gimbal Piston to stabilize relatively top-heavy payloads and quickly dissipate disturbances of the isolated table top.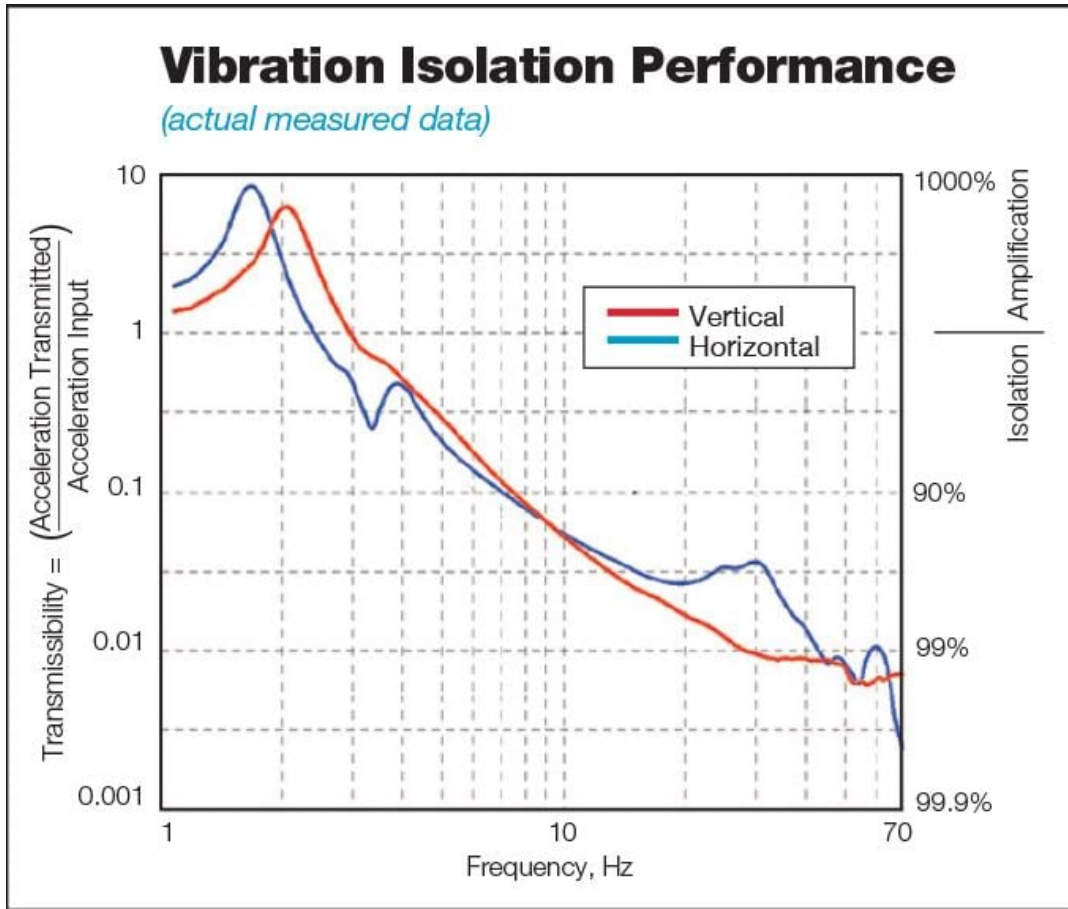

ERROR: There are no stock numbers to display!

Edmund Optics, Inc. - 101 E Gloucester Pike, Barrington, NJ 08007 | 800-363-1992 | Compliance@edmundoptics.com

Close Window



ERROR: There are no stock numbers to display!

Edmund Optics, Inc. - 101 E Gloucester Pike, Barrington, NJ 08007 | 800-363-1992 | Compliance@edmundoptics.com

Close Window

NOTES:

1. VIBRATION RATING:
5 Hz: Vertical = 70-85%, Horizontal = 75-90%
10 Hz: Vertical = 90-97%, Horizontal = 90-97%
2. FEED PRESSURE (psi): 80 (nitrogen or air)
3. CONSTRUCTION:
TOP: STAINLESS STEEL
CORE: ALL STEEL, HIGH DENSITY, CELL-SIZED HONEYCOMB CORE
FRAME: MEDIUM TEXTURE BLACK POWDER COAT
4. WEIGHT: 600 lbs
5. LOAD CAPACITY (lbs):
GROSS: 1400
NET: 350
6. ROHS: COMPLIANT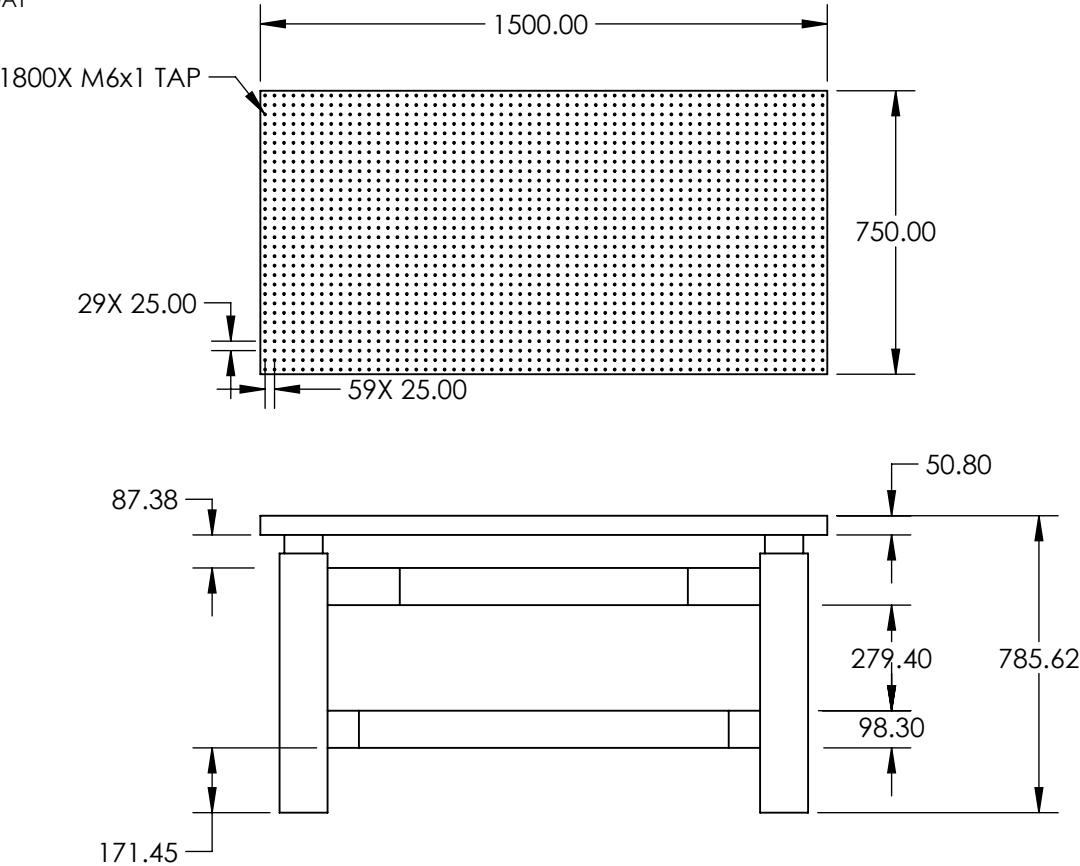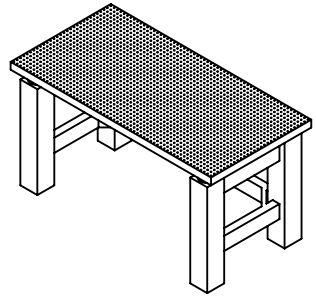

SPECIFICATIONS SUBJECT TO CHANGE WITHOUT NOTICE
DIMENSIONS ARE FOR REFERENCE ONLY

**FOR INFORMATION ONLY:
DO NOT MANUFACTURE
PARTS TO THIS DRAWING**

EO[®] Edmund Optics[®]

THIRD ANGLE PROJECTION		TITLE	750 x 1500mm, LOW PROFILE LAB TABLE	
ALL DIMS IN	mm	DWG NO	33782	SHEET 1 OF 1